295 DeKalb Pike, North Wales, PA 19454


TEL: 215-699-8700 FAX: 215-699-0370

Toll Free FAX: 1-800-255-4017

"L" SERIES MIXER

ASSEMBLY INSTRUCTIONS AND PARTS LIST

MIXER MODELS: L-1-300, L-1-CL, L-1-BK

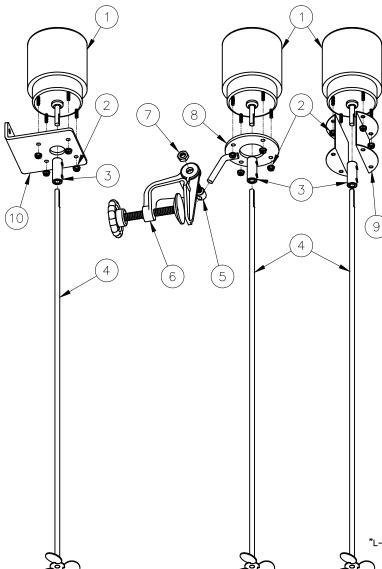

ATTENTION: TO INSURE SAFE AND EASY USE OF YOUR MIXER, READ THESE INSTRUCTIONS BEFORE USING.

Description

Laboratory mixers are designed for mixing of fluid in a container. The mixer is mounted using a supplied clamp or bracket, which is attached to a tank. The shafts with propeller are removable for cleaning.

General Safety Information

Warning: These mixers are intended for specific functions.

Please follow all instructions to insure
safe and proper use.

- (1) Be certain that all mounting hardware and the setscrews of coupling are properly tightened prior to operation.
- (2) Be certain propeller is not touching bottom or side of the mixing container before starting mixer.
- (3) Do not attempt to remove the mixing container or mixer when mixer is operating, as this could result in personal injury or damage to the mixer.
- (4) Always wear safety goggles when operating the mixer.

Assembly Instructions

- (1) Assemble Mixer as per the drawings. Make sure all hardware is tight.
- (2) Attach the coupling (Part # 3) to the air motor shaft (Part #1)
- (3) Insert the shaft and propeller assembly into the coupling. Tighten the setscrew.

"L-1-BK" MIXER

"L-1-300" MIXER MIXER MODELS
"L-1-CL" & "L-1-E"

| | | | MIXER MODELS | | | | | | |
|-------------|--|------|--------------------|-----------------------|-----------------------|-----------------------|-----------------------|--------------------|-------------------------|
| ITEM NO. | DESCRIPTION | QTY. | L-1-CL PART NO. | L-1-CL 24 PART NO. | L-1-CL 18 PART NO. | L-1-CL 12 PART NO. | L-1-E PART NO. | L-1-BK PART NO. | L-1-300 PART NO. |
| 1 | MIXER MOTOR | 1 | 100783 | 100783 | 100783 | 100783 | 100783 | 100783 | 100783 |
| 2 | #8-32 KEP NUT | 4 | 101221 | 101221 | 101221 | 101221 | 101221 | 101221 | 101221 |
| 3 | SHAFT COUPLING WITH SETSCREWS | 1 | 000967 | 000967 | 000967 | 000967 | 000967 | 000967 | 000967 |
| 4 | SHAFT ASS'Y. 5/16" DIA. X 30" LG. & 3" PROP. | 1 | 001073 | - | • | | - | 001073 | 001073 |
| | SHAFT ASS'Y. 5/16" DIA. X 24" LG. & 3" PROP. | 1 | • | 003129 | • | - | - | - | - |
| | SHAFT ASS'Y. 5/16" DIA. X 18" LG. & 3" PROP. | 1 | ı | - | 003128 | - | - | 1 | - |
| | SHAFT ASS'Y. 5/16" DIA. X 12" LG. & 3" PROP. | 1 | - | - | - | 004113 | - | - | - |
| | SHAFT ASS'Y. 5/16" DIA. X 30" LG. & 3" PROP. | 1 | - | - | - | - | 000970 | - | - |
| 5 | PIVOT BOLT | 1 | 000528 | 000528 | 000528 | 000528 | 000528 | - | - |
| 6 | CLAMP BODY ASSEMBLY | 1 | 002092 | 002092 | 002092 | 002092 | 002092 | - | - |
| 7 | 3/8-16 HEX NUT | 1 | 100642 | 100642 | 100642 | 100642 | 100642 | - | - |
| 8 | MOUNTING BRACKET ASSEMBLY | 1 | 000529 | 000529 | 000529 | 000529 | 000529 | - | - |
| 9 | MOUNTING BRACKET | 2 | - | - | - | - | - | 000968 | - |
| 10 | MOUNTING BRACKET | 1 | - | - | - | - | - | - | 003287 |